Singlet-Oxygen Measurements using the Fluorolog® System

Introduction

The chemistry, spectroscopy, and reactivity of singlet oxygen, $^{1}O_{2}$, is of considerable interest in environmental applications, novel organic syntheses and mechanisms of photo-oxygenation. Recently, research has been directed towards understanding the biological effects of $^{1}O_{2}$, such as the photodynamic destruction of cells in the presence of a dye, oxygen, and light.

understanding mechanisms that produce ¹O₂ is needed for the synthesis of new anti-cancer agents used in photodynamic therapy. Spectroscopic properties of the dye and characterized may be using fluorescence spectroscopy. The near-IR emission of ¹O₂ is around 1280 nm. The Fluorolog® system easily detects this appropriate signal with emission gratings and suitable detector. The standard R928P photomultipler tube responds to photons up to 800 nm. InGaAs detectors extend this range to nearly 1700 nm.

The DSS-IGA020L InGaAs detector is an excellent choice for $^{1}O_{2}$ measurements. It operates at liquid- N_{2} temperature to minimize thermal noise. Usable from 800–1700 nm, the InGaAs detector provides a smooth overlap with the R928P photomultiplier tube.

Experimental method and results

To demonstrate ${}^{1}O_{2}$ detection generated using a transition-metal complex, tris-2'-2'-bipyridine ruthen-ium(II) chloride, [Ru(bpy)₃]Cl₂, in D₂O,

was used as a $^{1}O_{2}$ generator. 2 The optical density of the solution was 0.3 A over the $D_{2}O$ baseline at 450 nm. To enhance the $^{1}O_{2}$ yield, the solution was saturated with O_{2} by bubbling pure O_{2} through it for 5 min before measurement. The λ_{max} of the $[Ru(bpy)_{3}]^{2+}$ complex is near 450 nm.

A Fluorolog[®] spectrofluorometer, equipped with a 450-W xenon lamp, single-grating monochromators. near-IR accessories (DSS-IGA020L InGaAs detector and emission grating at 600 grooves mm⁻¹, blaze = 1000 nm) was used to measure the ¹O₂ emission spectrum. A high-pass filter (λ = 780 nm) acted as an order-sorter, necessary when such large wavelength-ranges are spanned. Fig. 1 shows the emission spectrum with an integration time of 0.1 s and a bandpass of 12 nm. Note the excellent signal-to-noise ratio.

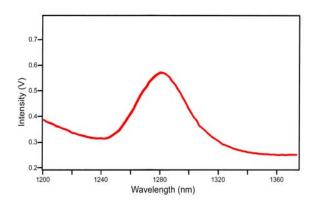


Fig. 1. Emission spectrum of ${}^{1}O_{2}$ generated by $[Ru(bpy)_{3}]Cl_{2}$ in $D_{2}O$.

Copyright © 2007 HORIBA Jobin Yvon; version 1.0

HORIBA

¹ Wasserman, H.H.; Murray, R.W., in *Singlet Oxygen*; Academic Press: New York, 1979.

² Mulazzani, *et. al. J. Phys. Chem.* **1994**, 98, 1145.



HORIBAJOBIN YVON

USA:

France:

HORIBA Jobin Yvon Inc., 3880 Park Avenue, Edison, NJ 08820-3012, Toll-Free: +1-866-jobinyvon Tel: +1-732-494-8660, Fax: +1-732-549-5125, E-mail: info@jobinyvon.com, www.jobinyvon.com HORIBA Jobin Yvon S.A.S., 16-18, rue du Canal, 91165 Longjumeau Cedex, Tel: +33 (0) 1 64 54 13 00, Fax: +33 (0) 1 69 09 93 19, www.jobinyvon.fr HORIBA Ltd., JY Optical Sales Dept, Higashi-Kanda, Daiji Building, 1-7-8 Higashi-Kanda Chiyoda-ku, Tokyo 101-0031, Tel: +81 (0) 3 3861 8231, www.jyhoriba.jp +49 (0) 89 462317-0 | Italy: +39 0 2 57603050 | UK: +44 (0) 20 8204 8142 Japan: Germany:

China: