


Symphony

1024 x 256 Thermoelectric Back-Illuminated Deep Depletion CCD Detector

The Most Sensitive Detector in the NIR with Minimized Etaloning Effects

The Jobin Yvon Back Illuminated Deep Depletion 1024 x 256 CCD is the best choice for low light level applications in the NIR (500 nm –1000 nm) including Raman, Photoluminescence, and Fluorescence spectroscopy. It offers the advantages of the high quantum efficiency from back illumination and the Deep Depletion technology for minimizing etaloning effects (oscillations superimposed on the CCD data due to the layer thickness of the classical Back Illuminated chip)

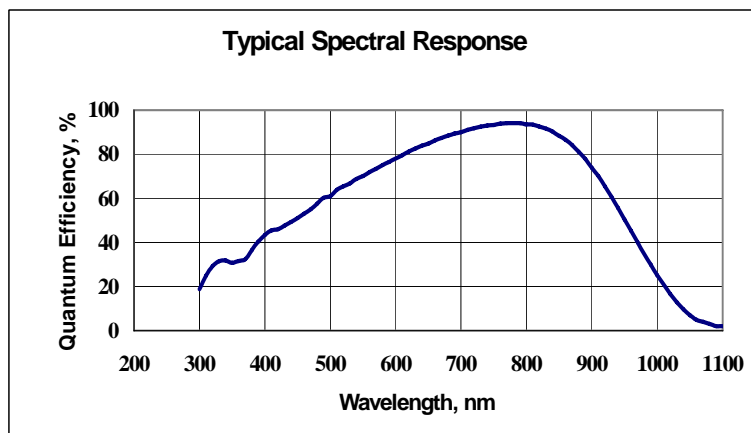


Features	Benefits
Scientific Grade 1 CCD	Ideally suited for low light level detection in a variety of spectroscopic applications
Back Illuminated Deep Depletion Technology	Enhanced NIR response and reduced etaloning
Liquid Nitrogen Cooling	Extremely low dark signal operation for extended integration times required with low signal levels
Excellent Linearity	Increased accuracy of data over the full dynamic range
Software Selectable Scan Rates	Optimize an experiment for the best combination of speed and sensitivity
Ethernet Connection to Host PC	Standard, easy to use interface with 100% data integrity
HORIBA Jobin Yvon's SynerJY™ Software	Complete control of a Symphony CCD and HORIBA Jobin Yvon Spectrograph system with full analysis capabilities
LabVIEW VIs and SDK Available	Flexible software to integrate a Symphony CCD into existing apparatus or as an OEM component

Specifications				
CCD Format		1024 x 256, Back Illuminated, Scientific Grade 1		
Pixel Size		26 μm x 26 μm		
Image Area		26.6 mm x 6.7 mm, 100% Fill Factor		
Cooling System		4 Stage Thermoelectric Cooling		
		Minimum	Typical	Maximum
Readout Noise	20 kHz		3.4 e ⁻ rms	6 e ⁻ rms
	1 MHz		20 e ⁻ rms	25x e ⁻ rms
Pixel Well Capacity		400 ke ⁻	700 ke ⁻	
Register Well Capacity			1000 ke ⁻	
Dark Current			2 e ⁻ /pixel/s	3 e ⁻ /pixel/s
Nonlinearity		< 0.4 % at 20 kHz scan rate < 1 % at all other scan rates		
Scan Rates		Software Selectable from 20 kHz to 1 MHz		
Software Selectable Gains		5 Software Selectable Gains		
Dynamic Range		16 bits		
Vertical Shift Rate		48 μs, 24 μs, 8 μs ¹		
Maximum Spectral Rate	20 kHz	14 Hz		
	1 MHz	182 Hz ^{1,2}		

Specifications subject to change without notice.

Mechanical Dimensions

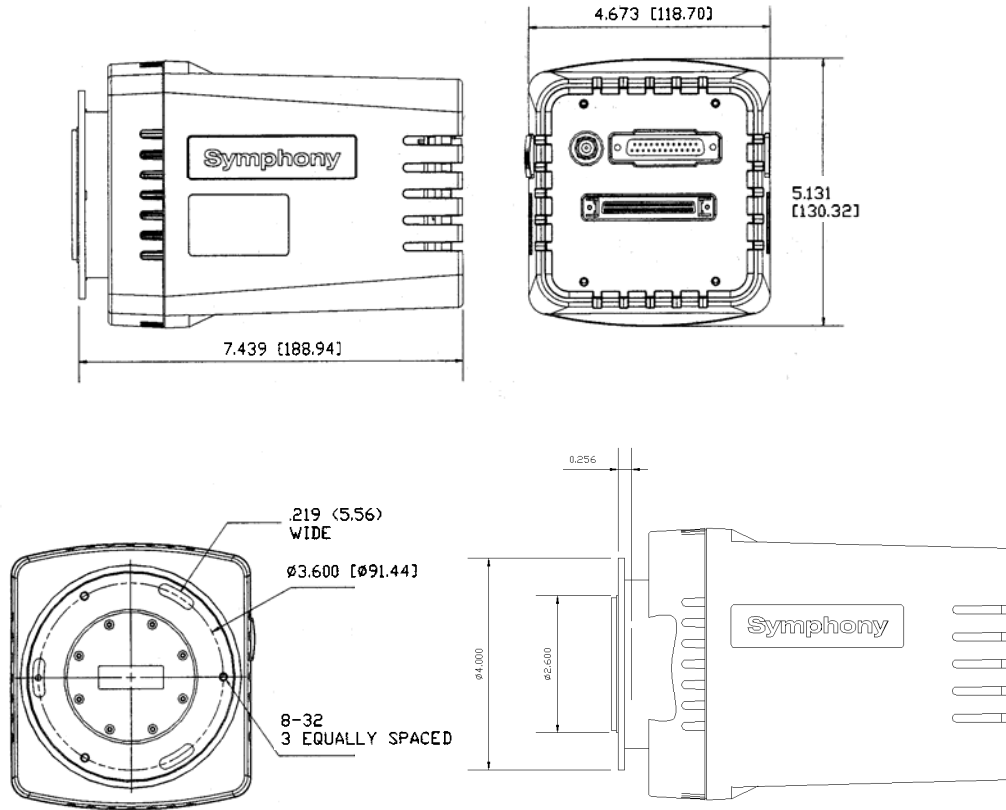

Units: Inches (mm)

Ordering Information:

CCD-1024x256-BIDD-STE Super Thermoelectric Cooled CCD System

Notes:

- 1 CCDs are guaranteed to have full Charge Transfer Efficiency at our standard shift rate of 48 μ s. At faster shift rates, a decrease in CTE may be observed
- 2 Highest Spectral rates are achieved when using the 1MHz ADC and a Parallel Transfer Time of 8 μ s